

Instructions to setup SoftLinX for Remote Access on Windows XP

Turn off Windows Firewall on both computers:

1. Click Start. Click Control Panel
2. If Control Panel is in Category View Click Network & Internet Connections
3. Click Network Connections
4. Right-click Local Area Network and select Properties
5. On the Advanced tab click Windows Firewall
6. Select Off
7. Click OK
8. Click OK

On the computer that will run the remote interface:

1. Click Start. Click Run... Type "dcomcnfg". Click OK
2. Expand Component Services, Computers, My Computer
3. Right-click My Computer and select Properties
4. Switch to the Default Properties tab
5. Check the checkbox label Enable Distributed COM on this computer
6. Under Default Authentication Level select Connect
7. Under Default Impersonation Level select Identify
8. Switch to the COM Security tab
9. Under Access Permissions click Edit Limits
10. Click Add
11. Type INTERACTIVE
12. Click OK
13. Check the Allow checkbox for all permissions
14. Click OK
15. Under Access Permissions click Edit Defaults
16. Repeat steps 10-14
17. Under Launch and Activation Permissions click Edit Limits
18. Repeat steps 10-14
19. Under Launch and Activation Permissions click Edit Defaults
20. Repeat steps 10-14
21. Click OK
22. Expand DCOM Config
23. Right-click SoftLinX.SLInterface and select Properties
24. On the Identity tab, select The Interactive User
25. Click OK
26. Restart the computer